

Investigation of Carryover or Cross-contamination in the Dionex ASE[®] 200

INTRODUCTION

Accelerated Solvent Extraction (ASE) is an innovative sample preparation technique that combines elevated temperatures and pressures with liquid solvents to achieve fast and efficient removal of analytes of interest from various matrices. With ASE, extractions can be done in very short periods of time and with minimal amounts of solvent as compared to conventional sample extraction techniques like Soxhlet or sonication. For example, 10-g samples can be completely extracted in less than 15 minutes with less than 15 mL of solvent. ASE has been demonstrated to be equivalent to existing extraction methodologies such as Soxhlet and automated Soxhlet for most RCRA (Resource Conservation and Recovery Act) analytes from solid and semisolid samples. It meets the requirements of U.S. EPA Method 3545, Pressurized Fluid Extraction.

With the small amount of solvent used relative to the sample size, carryover or cross-contamination could be potential concerns with ASE and the Dionex ASE 200. Two sets of experiments were conducted to investigate these concerns. The experiments performed included the extraction of heavily loaded soil and sediment samples followed by extracting blank samples and the determination of target analytes in both extracts.

INSTRUMENTATION

Dionex ASE 200 Accelerated Solvent Extractor
 Dionex DX 500 HPLC system with AD20 (UV detector) and fluorescence detector
 Gas chromatograph (GC) with electron capture detector (ECD)

EXPERIMENTAL

In the first experiment, 7-g samples of a highly contaminated soil (SRS100-103, Fisher Scientific, 11 wt% solvent extractable) were extracted at 17.2 MPa (2500 psi), 100 °C with toluene/methanol (1:1), with 5-minute heat-up and 5-minute static times. Analyses of the extracts were done by HPLC using UV at 254 nm and fluorescence at 325-nm excitation and 410-nm emission. In this case, the target analytes were polycyclic aromatic hydrocarbons (PAH). After the extraction, the cell was removed and replaced with a cell containing clean sand. A second extraction was performed using the same conditions. These steps were repeated with three separate new samples to determine reproducibility. No rinsing of the instrument was performed between the extractions.

The second experiment was conducted using Standard Reference Material (SRM) 1939 (river sediment from the National Institute of Standards and Technology (NIST), Gaithersburg, MD). This sample has a high level of coextractable components (about 30 wt%) in addition to the PCBs for which this material is certified. In this experiment, 8 g of the sediment were extracted with hexane/acetone (1:1) at 100 °C and 17.2 MPa (2500 psi) with 5-minute heat-up and 5-minute static periods.

Table 1 Carryover of PAH in the ASE 200 After Extraction From Contaminated Soil

	HPLC With UV Detection	HPLC With Fluorescence Detection	ASE Recovery vs. Certified Values
PAH Carryover n=3	0.024%	0.13%	106% Recovery 5.5% RSD

After the extraction, a second extraction was performed on the same sample and collected in a separate vial. This process was repeated with three samples. No rinsing of the instrument was performed between the extractions. The analysis of the extracts was performed by GC with ECD. These experiments were performed to determine not only if material remained in the tubing of the instrument, but to determine the completeness of the extraction of the sample in the cell itself.

RESULTS AND DISCUSSION

The experiments on the contaminated soil were performed to determine if any material remained and was not completely flushed from the tubing and valves that are in line with the extraction cell. Carryover was determined by comparing the total area counts from the samples to those obtained from the blanks for the UV and fluorescence detectors on the HPLC analysis method for PAH. The results are shown in Table 1. These are the percentages of the total area seen in the first extraction of the sample that were obtained from the extract of the empty cell. These results show that an insignificant level of carryover exists. It should be noted that it is unlikely that this carryover was due to analytes because the recovery of the PAH was quantitative compared to the certified values. No PAH were detected in second extracts or blanks (minimum detectable quantity (MDQ) was 0.2 mg/kg, and the certified concentration of the PAHs ranged from 31 to 1500 mg/kg).

In the experiments with the river sediment sample, the total area counts (minus the solvent blank) of the first extract were compared to those of the second. The results are summarized in Table 2. This small amount of material

Table 2 Carryover of PCBs in the ASE 200 After Extraction From Contaminated River Sediment

	GC Analysis With ECD	ASE Recovery vs. Certified Values
PCB Carryover n=3	2.6%	102% Recovery 4% RSD

in the second extract was actually residual coextractable material because the recovery of the target compounds, PCBs, was quantitative compared to the certified values. No PCBs were detected in the second extract (MDQ = 0.1 µg/kg, and the concentration of the analytes ranged from 180 to 3700 µg/kg). This indicates that the lines are cleaned out well, and complete extraction is achieved with these small volumes of solvent.

The ASE 200 has low dead-volume, which translates into no cross-contamination between samples. It can be programmed to perform automatic rinses between samples to allay concerns of cross contamination between samples. Extensive testing for this was also conducted as part of the EPA equivalency studies. ASE is an exhaustive extraction technique, and when quantitative extraction occurs, no analytes are left in the sample for subsequent extractions.

ASE is a trademark of Dionex Corporation.



Printed on recycled and recyclable paper with soy-based ink.

Dionex Corporation
1228 Titan Way
P.O. Box 3603
Sunnyvale, CA
94088-3603
(408) 737-0700

Dionex Corporation
Salt Lake City Technical Center
1515 West 2200 South, Suite A
Salt Lake City, UT
84119-1484
(801) 972-9292

Dionex U.S. Regional Offices
Sunnyvale, CA (408) 737-8522
Westmont, IL (630) 789-3660
Houston, TX (281) 847-5652
Atlanta, GA (770) 432-8100
Marlton, NJ (856) 596-0600

Dionex International Subsidiaries
Austria (01) 616 51 25 Belgium (015) 203800 Canada (905) 844-9650 France 01 39 46 08 40 Germany 06126-991-0
Italy (06) 66 51 50 52 Japan (06) 6885-1213 The Netherlands (0161) 43 43 03 Switzerland (062) 205 99 66 United Kingdom (01276) 691722
* Designed, developed, and manufactured under an NSAI registered ISO 9001 Quality System.
<http://www.dionex.com>



LPN 0766 5 M 7/96
© 1996 Dionex Corporation